RAPTOR

SERVER ANALYSIS SOFTWARE

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RAPTOR

SERVER ANALYSIS SOFTWARE

Welcome to RAPTOR SERVER Analysis Software, the analysis software for the [ALEX multiplex test](#). Please log in below to continue.

<p>Log in with Azure AD / Office 365</p> <p>If your organization integrates the software into its Microsoft Azure/Office 365 user directory (single-sign-on) click the button below. You will be redirected to the Microsoft login screen.</p> <p> Log in</p>	<p>Log in with Username & password</p> <p>Please enter your username and password.</p> <p>Your username</p> <p>Your password </p> <p> Log in</p> <p>Forgot your password? No problem, you can request a new one.</p>
----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------	-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------

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Log in with Username & password

Please enter your username and password.

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Contract approval required

Please review the following documents. You are required to confirm that you have read and agree with each document in order to use RAPTOR SERVER:

- [Bitte lesen](#) (last updated 23.11.2023 18:00)
- [RAPTOR SERVER IFU](#) (last updated 04.12.2023 11:05)
- [Terms of Service](#) (last updated 22.11.2023 01:00)

Confirm

By clicking the button, I confirm that I have reviewed these documents and agree.

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QC	Product type / Source	QR-Code	Sample Code	Analysis Date	
❌ Failed	Research MAX 45k	99AFG0F7	NK FOX	30/03/2023 20:50	Details Report
❌ Failed	Research MAX 45k	99AFG0F8	N.G.	30/03/2023 20:50	Details Report
❌ Failed	Research MAX 45k	99AFG0F9	E18735	30/03/2023 20:50	Details Report
❌ Failed	Research MAX 45k	99AFG0FA	20220325w56	30/03/2023 20:50	Details Report
❌ Failed	Research MAX 45k	99AFG0EE	HM100038	30/03/2023 20:50	Details Report
❌ Failed	Research MAX 45k	99AFG04D	786118	30/03/2023 20:37	Details Report
❌ Failed	Research MAX 45k	99AFG016	HM100225	30/03/2023 20:37	Details Report
❌ Failed	Research MAX 45k	99AFG0CA	E17685_M	30/03/2023 20:37	Details Report
❌ Failed	Research MAX 45k	99AFG0CC	E17685_M	30/03/2023 20:37	Details Report
❌ Failed	Research MAX 45k	99AFG015	E17685_F	30/03/2023 20:37	Details Report

In total there are 270 new measurements in the selected time period.

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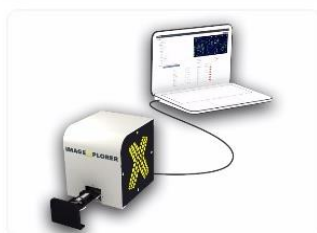
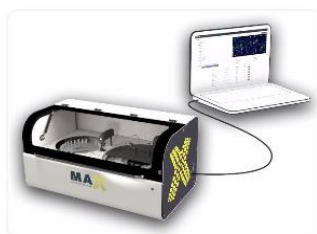
9K-Test-Instrument

Not initialized

[Initialize](#) [Shutdown](#) [Monthly Maintenance](#)

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Name:

E-Mail address:

Language: English (en)
The language is used to determine the display language.

Culture: English (United States) (en)
The culture is used to format numbers and dates.

Timezone: (UTC+01:00) Amsterdam, Berlin, Bern, Rome, Stockholm, Vienna
Configure your timezone to see date and times in your local time instead of UTC.

Multi-Factor Authentication

[Manage MFA](#)


Use multi-factor authentication (MFA) to significantly increase security. It is highly recommended to enable MFA; this adds an additional layer of protection against phishing attacks or weak passwords.

Configuration

Research spots:
If enabled, "research use only" arrays. This does not affect the

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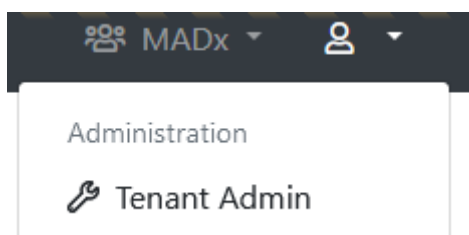General configuration

*** Name:** MADx

Default timezone: (UTC+01:00) Amsterdam, Berlin, Bern, Rome ▾
Configure a timezone to see date and times in (instead of UTC). Please note that users can choose to use a different timezone.

Default report template (ALEX2): ALEX2 ▾

Default report template (ALEX3, ALEX Air, ALEX Food): ALEX (modern) ▾

Default report template (FOX): FOX (legacy) ▾

Default report template (PAX): PAX ▾

Default report language: English (en) ▾

Default ALEX2 MOD: < No MOD > ▾

Default FOX MOD: < No MOD > ▾

Default PAX MOD: < No MOD > ▾

Force MOD: False ▾
Requires a measurement to always have an MOD assigned. If this setting is enabled, a measurement cannot be approved without an MOD assigned.

Hide Global MOD: False ▾
If this setting is enabled, only the tenant-specific MOD will be available to select, and the global system-wide MOD will be hidden.

Display MOD Excluded Spots As Research Use Only: False ▾
If this setting is enabled, spots excluded by an MOD will not be hidden from the measurement image, but displayed as a "RUO spot".

MAX device default settings

*** Default Minimum RunId:** 10 ▾
When a new MAX device is created, this value will be used for its **Minimum Run Id**

Tenant-specific batch releases

Restrict this tenant to only allow the following batches to be scanned or to restrict certain batches. Enter one per line.

Use:

Whitelist

Blacklist

Whitelisted Batches:
Enter one batch code including product identifier per line, for example "02AAA"

Blacklisted Batches:
Enter one batch code including product identifier per line, for example "02AAA"

Company Details

*** Legal Entity Name:** MacroArray Diagnostics

*** Street and Number:** Street 4

*** Zip Code:** 1230

*** City:** New York

*** Country:** United States of America ▾

*** Phone:** +4318652573

*** E-Mail:** office@macroarraydx.com

*** Legal Representative:** Dr. Christian Harwanegg

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Tenant-specific batch releases

Restrict this tenant to only allow the following batches to be scanned or to restrict certain batches. Enter one per line.

Use:

Whitelist

Blacklist

Whitelisted Batches:

02AAA
02ABC
02BAA

Enter one batch code including product identifier per line, for example "02AAA"

Blacklisted Batches:

Enter one batch code including product identifier per line, for example "02AAA"

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You requested a new measurement from ImageExplorer "IX Trainingsraum". Please wait until the measurement is completed ...

Invalid batch detected! The measurement can't be processed because its batch is blocked.

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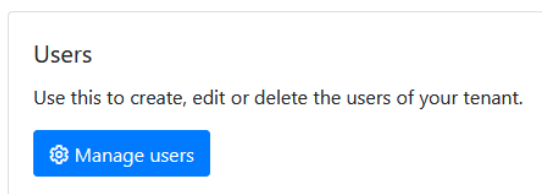
Login Credentials

Type:

Please select if the user should be validated against Azure AD (Office 365), or if you want to specify username and password yourself. It is recommended to use Azure AD whenever possible. This allows for increased security by supporting two-factor authentication. Also, it is more convenient for the user, because he or she does not have to remember a new username and password (single-sign-on).

Username:

It is a good idea to use the users e-mail-address as his username, because the username needs to be unique across the entire system.

Password:

Access Rights

Preconfigured groups

Use these pre-configured groups to automatically assign a set of access rights to this user. You can still add or remove specific access rights.

[Reset](#) [Lab User](#) [Super Lab User](#) [Support/Distributor User](#) [System Administrator](#)

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Tenants

This user is connected to the following tenants and has the following access rights:

<input type="checkbox"/> Tenants	Can share measurements	Can edit measurements	Can edit personal data	Can approve measurements	Can manage basic tenant configuration	Can manage API keys	Can manage ImageXplorers	Can view agent connections	Can manage users	Can manage users of Result Sharing Portal	Can manage report templates	Can manage measurement data completion	System Administration	Can unapprove measurements	Can use a custom calibration string	Can manage Run Configurati	Can skip r
<input type="checkbox"/> Another MADx Tenant	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>
<input type="checkbox"/> MADx Tenant	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>

Additionally this user can be added to the following tenants by checking the tenant name and hitting save afterwards. User rights w applied from current open tenant user.

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Additionally this user can be added to the following tenants by checking the tenant name and hitting save afterwards. User right applied from current open tenant user.

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Tenants

This user is connected to the following tenants and has the following access rights:

[Edit selected tenants rights](#)

<input type="checkbox"/>	Tenants	Can share measurements	Can edit measurements	Can edit personal data	Can approve measurements	Can manage basic tenant configuration	Can manage API keys	Can manage Imagey
<input checked="" type="checkbox"/>	Another MADx Tenant	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>

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General information

Name:

Key:
This key is required when installing an Agent.

ImageExplorer Serialnumber:
Serialnumber of connected ImageExplorer.

Notes:

Internal Notes:
For internal use only.

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Analysis Log

```
'PreProcessedPixels' took 128ms.
Using definition '02_ALEXII_IVD_v1.6.0'.
'ApplyArrayDefinition' took 74ms.
Parsing calibration string ...
'ApplyDefinitions' took 75ms.
'FindGrid' took 1ms.
'AlignFeatures' took 862ms.
'CropImage' took 1ms.
'QuantifyFeatures' took 233ms.
'SubtractBackground' took 681ms.
Parsing calibration string ...
Using calibrators from QR-Code.
Total IgE (tIgE) normalized by calculating 0,330999755859375E-07x*+0,05697999954223633x+-30,69269921875
(x=8515), which results in 515.
'ReQuantify' took 57ms.
```

QC Log

These are the values used for QC (ALEX v2): Curve Correction=0,9997370867894242, tIge=8515, GD1=28985 (SNR=185.38), GD2=29369 (SNR=185.28), GD3=28569 (SNR=147.53), GD4=28305 (SNR=128.1)
Reference values are as follows: Guide Dots must have a value above 5000; each Guide Dot spot must exist and be unique; Guide Dot spot must not be set to bad. Average background must be between 4000 and 16000.
QC (ALEX v2) succeeded.

Background Log

Calculated background statistics for 164 pixels (from a total of 182, skipped smallest 9, and highest 9 pixels). The average is 9.419, the minimum 7973 and the maximum 11792 (minimum and maximum is calculated including skipped pixels). This results in an average difference of 628, which is an SD of 49,17 and a CV of 0,52%.
For comparison, there is an overall total of 282944 pixels. The overall average is 14,071, the minimum 6,012 and the maximum 64,965.

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New Measurements

Search by QR-Code, patient, sample, tag, analysis date, approval date ...

Selected measurements

<input type="checkbox"/>	Source	QC	Tags	Product type
<input type="checkbox"/>	IX	✓ OK		ALEX2
<input checked="" type="checkbox"/>	IX	✓ OK		ALEX2
<input checked="" type="checkbox"/>	IX	✓ OK		ALEX2
<input checked="" type="checkbox"/>	IX	✓ OK		ALEX2

- Download CSV format
- Download Excel format
- Approve
- Update interpretation
- Choose Reporttemplate
- Modify Tags**
- Move
- Delete

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Re-process measurements

You are about to re-process the following measurements. undone, the existing analysis result will be irreversibly over

QR-Code	Sample Code
02APK3C8	Dummy8
02APK258	Dummy1
02APK07F	Dummy13

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New Measurements

Search by QR-Code, patient, sample, tag, analysis date, approval date ...

<input type="checkbox"/>	Source	QC	Tags	Product type	QR-Code	Sample Code
<input type="checkbox"/>	IX	✓ OK		ALEX2	02AGT957	
<input type="checkbox"/>	IV	✓ OK				

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New Measurements

Search by QR-Code, patient, sample, tag, analysis date, approval date ...

<input type="checkbox"/>	Source	QC	Tags	Product type
<input type="checkbox"/>	Recaptured	✓ OK	recaptured	ALEX1
<input type="checkbox"/>	⚠ Recaptured	✓ OK	recaptured	ALEX1
<input type="checkbox"/>	QR-Code was scanned more than once.		recaptured	ALEX1

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RAPTOR New measurements Approved measur

New Measurements

Search by QR-Code, patient, sample, tag, analysis date, approval date ...

Selected measurements

<input type="checkbox"/>	Source	QC	Tags	Product type	QR-Code
<input checked="" type="checkbox"/>	IX	✓ OK		ALEX2	02AMC180
<input checked="" type="checkbox"/>	IX	✓ OK		ALEX2	02AMC078
<input checked="" type="checkbox"/>	IX	✓ OK		ALEX2	02AKV2D9

- Download CSV format
- Download Excel format
- Approve**
- Update interpretation
- Re-Process
- Move
- Delete

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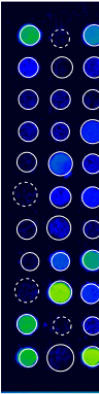
Approve measurements

You are about to approve the following measurements.

There are measurements with one or more warnings. Please make sure to double-check these measurements before approval.

QR-Code	Sample Code	Analysis Date
02AKV035		28.04.2022 11:14

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Export requested	# of measurements included	Export completed
8/16/2021 11:07 AM	4	0 / 4
8/16/2021 11:06 AM	4	8/16/2021 11:06 AM
8/16/2021 11:04 AM	3	8/16/2021 11:04 AM

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MADx Tenant

- Administration
 - Tenant Admin
 - System Admin
 - Exports**
 - Personal Settings
 - System Notifications
 - Log out
- Support
- Imprint

14.02.2024 16:15

Exports

Certain types of data exports (for example, the export of many PDF reports at a time) are possible. Please note that you can keep working, all exports are done in the background.

Export multiple measurements

Export requested	# of measurements included
06.07.2022 08:52	3

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Export multiple measurements

Please note that **Report** and **ASTM** export is only available for certain product types.

Product Type
ALEX2

Analysis Date From
TT . MM . JJJJ

Tag
use , to separate tags

Export CSV

QR Code

Analysis Date
TT . MM . J

Approved or not
New meas

Export Excel

Search Measurements **Cancel**

Export multiple measurements

Please note that **Report** and **ASTM** export is only available for certain product types.

Product Type

ALEX2

QR Code

Analysis Date From

TT . MM . JJJJ

Analysis Date To

TT . MM . JJJJ

Tag

Approved by

Approved

use , to separate tags

Export CSV

Export

 Start export ...

← Back

122 Measurements found that match your criteria

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Configure report template "New report template"

The screenshot shows the configuration interface for a new report template. It has four tabs: General, Design, Header, and Content. The 'General' tab is active. It contains three main settings:

- Name:** A text input field containing "New report template". A tooltip explains: "This is the name of the template which shows up in the dropdown list while generating a report."
- Base Template:** A dropdown menu showing "Version 2 (ALEX2 only)". A tooltip explains: "A base template which specifies basic functionality must be chosen."
- Preferred language:** A dropdown menu showing "English (United States) (en-us)". A tooltip explains: "A default language can be specified which will be used if there isn't yet a language stored to a measurement's report."

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PDF naming

Setting Type: Manual

The filename of the report PDF can be customized. Either built automatically based on your selection below or manually adjusted in the order of your needs.

Name Pattern: Report_{PatientId}_{MeasurementQrCode}

Available Elements: {PatientId} ; {PatientName} ; {SampleId} ; {MeasurementQrCode} ; {DateAndTime}

PDF naming

Setting Type: Automatic

The filename of the report PDF can be customized. Either built automatically based on your selection below or manually adjusted in the order of your needs.

Include Patient ID:

Include Patient Name:

Include Sample ID:

Include Measurement QR-Code:

Include Date & Time:

Footer

Footer text:

Will be printed in the footer if provided and rendered as HTML.

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Configure report template "New report template"

The screenshot shows the 'Design' tab of the 'Configure report template' interface. It contains three main settings:

- PDF format:** Radio buttons for 'A4' (selected) and 'Letter'. A note states: 'Specifies the paper size of the output PDF.'
- Font:** A dropdown menu showing 'Arial'. A note states: 'The font family of the whole report texts can be adjusted.'
- Background image:** A text input field followed by a 'Browse' button. A note states: 'If set, this image is used as background on the first page of the PDF.' Below this field, it says 'Maximum file size 1 MB'.

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Logo (left):

Maximum file size 1 MB

This logo will appear on the **top left corner** of the first page of the report.

Text (center):

This text will appear either on the left, right, or center of the top of the first page (depending on your configured logos). Supports the placeholders `{{ PATIENT_PORTAL_QRCODE }}` and `{{ PATIENT_PORTAL_URL }}` to dynamically insert measurement information.

Logo (right):

Maximum file size 1 MB

This logo will appear on the **top right corner** of the first page of the report.

Small logo:

This logo is used on all detail pages. If not set, the default logo is used.

Signatures

Left:

Maximum file size 1 MB

This signature or image will appear on the last page, left aligned.

Center:

Maximum file size 1 MB

This signature or image will appear on the last page, center aligned.

Right:

Maximum file size 1 MB

This signature or image will appear on the last page, right aligned.

Indicators

Custom intensity indicator:

A custom SVG for displaying allergen intensity can be pasted here. Will be styled according to specified colors.

The SVG element must contain the attributes **width** and **height** with a minimum dimension of 16 for correct display in PDF.

SVGs paths which should be colored must require the **colorize** class as well as a fill defined in the **defs** segment of the SVG.

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Configure report template "New report template"

Header content:		Left half	Right half
		Custom headline	Custom headline
	<input type="checkbox"/>	Patient Name	Doctor Information
	<input checked="" type="checkbox"/>	Patient Name	None
	<input type="checkbox"/>	Date of Birth	None
	<input type="checkbox"/>	Sample ID	None
	<input type="checkbox"/>	QR-Code	Additional Info text
	<input type="checkbox"/>	Analysis Date	None
	<input type="checkbox"/>	Amount of tested allergens	None
	<input type="checkbox"/>	Test method	Quality Control text

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General	Design	Header	Content
Allergens to include:	All allergens		Specifies which kind of allergens should be included in the report.
Allergens sort order:	Alphabetically by allergen		Specifies the sort order of the allergens.
Include Overview:	<input checked="" type="checkbox"/>		Whether the allergen overview should be printed or not.
Include product type in overview:	<input checked="" type="checkbox"/>		Whether the product type should be printed in the header of allergen overview.
Include Total IgE:	On the front page		Whether the information of total IgE should be included in the report and where.
Cross-reactive information:	Include text		This section will be displayed underneath the allergen list itself.
Include Interpretation:	Include manually		It can be specified if the Raven report interpretation information should be included in the report or not. This section is printed underneath the allergens overview.

i Additional information (below the list of results)	Those informations are integrated into the report in an extra section underneath the allergen list.
Approval date: <input type="checkbox"/>	
Assay date: <input type="checkbox"/>	
Doctor info: <input type="checkbox"/>	The information about the doctor which accompanied the measurement/patient.
Print date: <input checked="" type="checkbox"/>	When the report was generated.
Report info: <input type="checkbox"/>	The „additional info“ field belonging to the measurement.
Sample date: <input checked="" type="checkbox"/>	

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MADx Administration Tenant Admin

Tenant configuration

The general configuration settings for your tenant.

[Configure Tenant](#)

General configuration

*** Name:** MADx

Default timezone: (UTC+01:00) Amsterdam, Berlin, Bern, Rome, Stockholm, Vienna
Configure a timezone to see date and times in (instead of UTC). Please note that users can choose to use a different timezone.

Default report template (ALEX 1): Version 1 (ALEX1 & ALEX2)

Default report template (ALEX 2): Version 2 (ALEX2 only)

Default report language: English (United Kingdom) (en-gb)

Default MOD: < No MOD >

Require ImageXplorer testarray scan: Warn after 60 days, enforce after 90 days.

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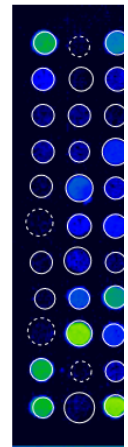
The screenshot shows the RAPTOR web interface. At the top, there are navigation tabs for 'New measurements' and 'Approved measurements'. Below this is a search bar and a 'Download selected measurements' dropdown menu. The main content area displays a table of 'Approved Measurements' with columns for Source, Tags, Product type, QR-Code, and Sample C. Three rows are visible, all with 'IX' as the source and 'ALEX2' as the product type. A dropdown menu is open over the table, showing options: 'Report (PDF format)', 'CSV format', 'Excel format', 'ASTM format', 'Update interpretation', 'Choose Reporttemplate', 'Share', 'Move', and 'Unapprove'. The 'Report (PDF format)' option is highlighted with a yellow circle.

	Source	Tags	Product type	QR-Code	Sample C
<input checked="" type="checkbox"/>	IX		ALEX2	02AKV39D	
<input checked="" type="checkbox"/>	IX		ALEX2	02AKV39D	
<input checked="" type="checkbox"/>	IX		ALEX2	02AKV39D	

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Tags:

MOD:

Report Template: Version 2 (ALEX2 only) - Deutsch (de)

Raptor [New measurements](#) [Approved measurements](#) [Start measurement](#) [Automats](#)

Report

Template:

Language:

[Save](#) [Use defaults](#)

Usually this tenants default template and language will be used when sharing or exporting measurements. You may change the template and language associated with this measurement by saving your current configuration.

[Download PDF](#)

[Manage report template and settings](#)

You have access to manage the report templates and settings via the tenant administration. Please note that any changes will affect all users.

Patient ID:	
Patient Name:	
Date of Birth:	
Sample ID:	
Barcode:	02AAE
Sampled on:	
Tested on:	12/18/
Printed on:	12/19/
Note:	The internal QC (Plausibility)

Lab report: Summary on de

Pollen	Grass Pollen
	Tree Pollen
	Weed Pollen
Mites	House Dust M
Microorganisms	Fungal Spores
	Legumes
	Grain
Plant-Based Food	Spices
	Fruits
	Vegetables
	Nuts & Seeds
	Milk

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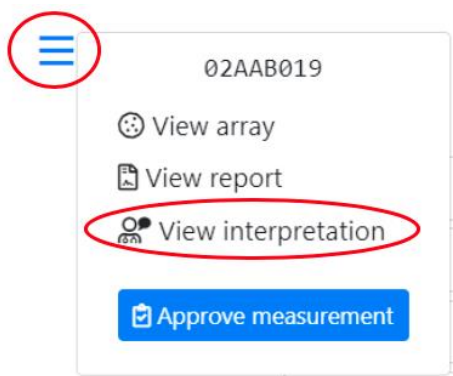
Raven Interpretation Summary

Sample Information

The sample was tested on ALEX[®] Barcode 02AAB019, interpretation date 12/19/2019.

Of the tested 295 allergens, 25 were above the detection limit of 0.3 kU_L/L. A sensitization can be an indicator of an allergy. For selected allergens, comments for interpretation are listed below.

Total IgE: 512 kU/L

The measured total IgE was 512 kU/L. Above a total IgE of 100 kU/L, allergy is likely.

Cross-Reactive Allergen Sensitization detected

Sensitization against allergen molecules which are markers of (broad) cross-reactivity between different allergen sources.

Detected cross-reactive allergen sensitizations:

- PR-10s: AIn g 1, Bet v 1, Cor a 1.0103, Fag s 1
- Poicacins: AIn g 4, Pnl p 7

PR-10 Proteins (PR10)

PR-10 inhalative: The major birch pollen allergen, Bet v 1, represents the prototype of all PR-10 allergens and is the primary sensitizer in regions with birch pollen exposure. The presence of PR-10 allergens in Fagales tree pollen explains IgE cross-reactivity between pollen from hazel, alder, beech, oak and hornbeam. PR-10 nutritive: PR-10 allergens in raw fruits, nuts, vegetable and legumes can induce oral allergy syndrome and sometimes severe allergic reactions in sensitized individuals. PR-10 allergens are not stable to heat and digestion.

Poicacins (Ca++ binding proteins, PC)

Members of the PC family can cause inhalative symptoms. PC allergens can be found in pollen from grasses, trees and weeds. Inhalative symptoms manifest as allergic rhino-conjunctivitis and/or allergic asthma.

Request new interpretation

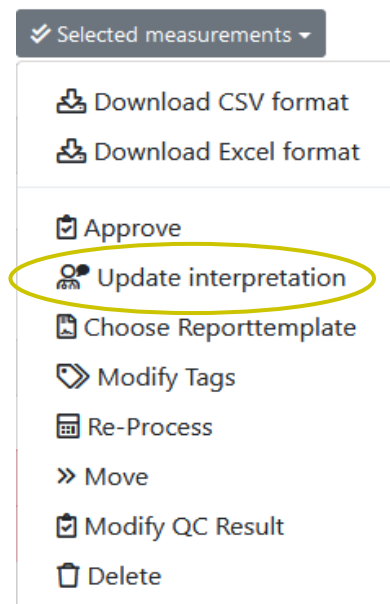
This measurement already has an interpretation stored. Usually, it should not be necessary to request those again, but it could be necessary when the patients information has changed.

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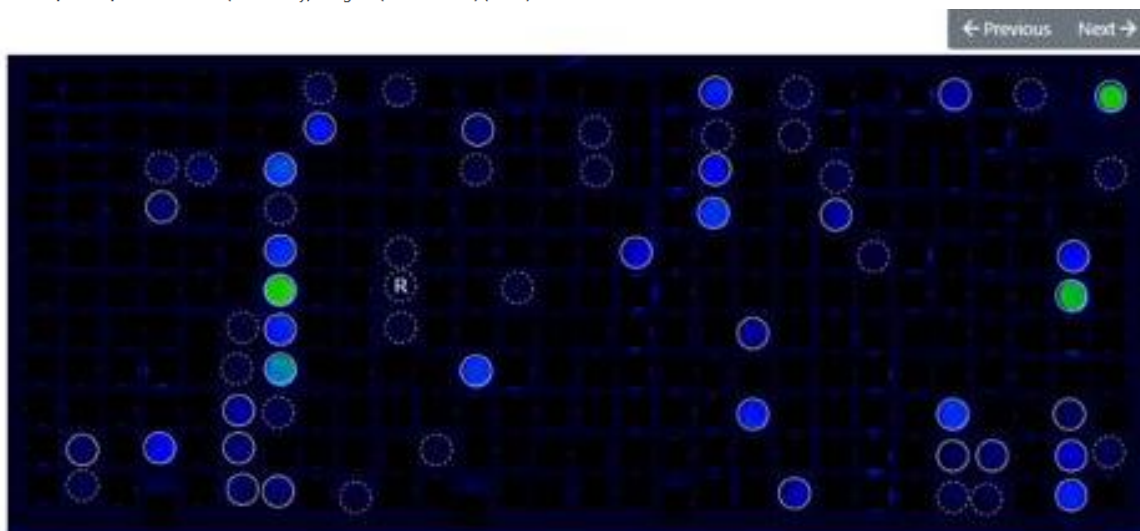
Tags:

MOD:

Report Template: Version 2 (ALEX2 only) - English (United States) (en-us)

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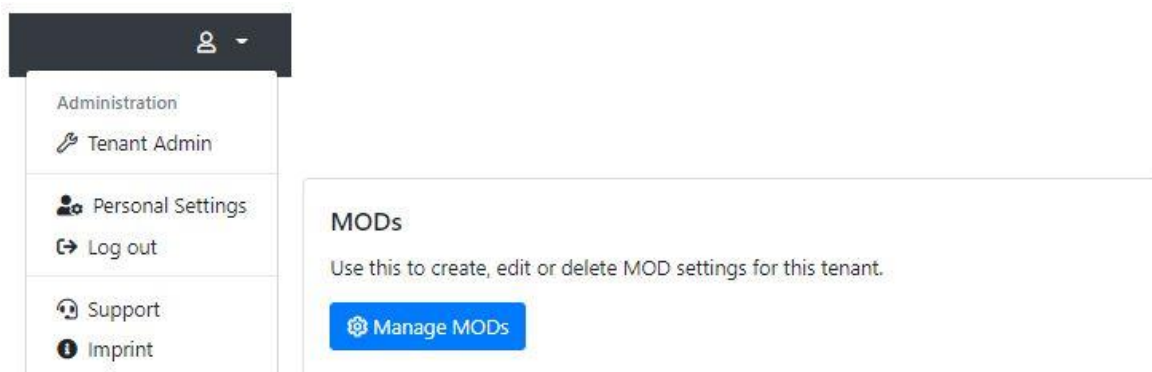
Raptor New measurements Approved measurements Start measurement Automats

Create new MOD

Name:

Type:

Deactivated:

You cannot delete an MOD once it is in use. But you can deactivate it, so that it's not available for new measurements anymore.

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* Name:

Default timezone:

Configure a timezone to see date and times in (instead of UTC). Please note that users can choose to use a different timezone.

Default report template (ALEX 1):

Default report template (ALEX 2):

Default report template (FOX):

Default report template (PAX):

Default report language:

Default ALEX2 MOD:

Default FOX MOD:

Default PAX MOD:

Force MOD:

Requires a measurement to always have an MOD assigned. If this setting is enabled, a measurement cannot be approved without an MOD assigned.

Hide Global MOD:

If this setting is enabled, only the tenant-specific MOD will be available to select, and the global system-wide MOD will be hidden.

Display MOD Excluded Spots As Research Use Only:

If this setting is enabled, spots excluded by an MOD will not be hidden from the measurement image, but displayed as a "RUO spot".

Require ImageXplorer ConfigXplorer scan: Warn after 60 days, enforce after 90 days.

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Global_MOD_ALEX_v1.17

MOD_ALEX_v1.17

Force MOD: True
Requires a measurement to always have an MOD assigned. If this setting is enabled, a measurement cannot be approved without an MOD assigned.

Hide Global MOD: True
If this setting is enabled, only the tenant-specific MOD will be available to select, and the global system-wide MOD will be hidden.

Display MOD Excluded Spots As Research Use Only: False
If this setting is enabled, spots excluded by an MOD will not be hidden from the measurement image, but displayed as a "RUO spot".

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
Force MOD:	False	◆
Requires a measurement to always have an MOD assigned. If this setting is enabled, a measurement cannot be approved without an MOD assigned.		
Hide Global MOD:	False	◆
If this setting is enabled, only the tenant-specific MOD will be available to select, and the global system-wide MOD will be hidden.		
Display MOD Excluded Spots As Research Use Only:	True	◆
If this setting is enabled, spots excluded by an MOD will not be hidden from the measurement image, but displayed as a "RUO spot".		

Configuration

Calibration string:	<input type="text"/>
Allows to override the calibration information from the cartridge's QR-Code. Leave empty to use the calibration information from the QR-Code on the cartridge (the usual behaviour).	
Research spots:	Show research spots ◆
If enabled, "research use only" spots will be displayed on measurement details and arrays. This does not affect the export behaviour or the reports of measurements.	

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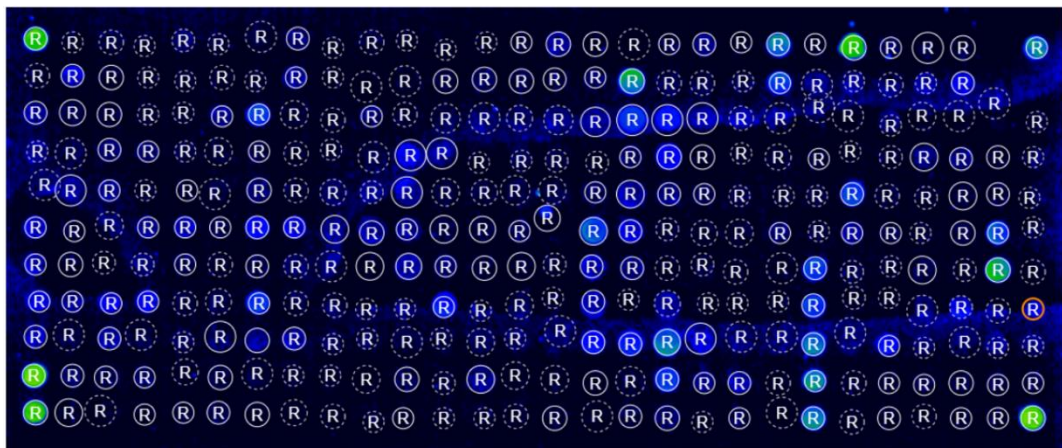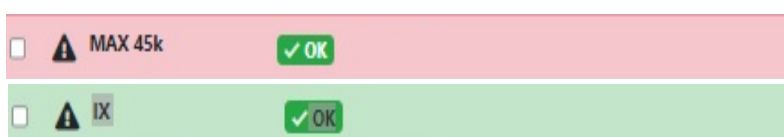
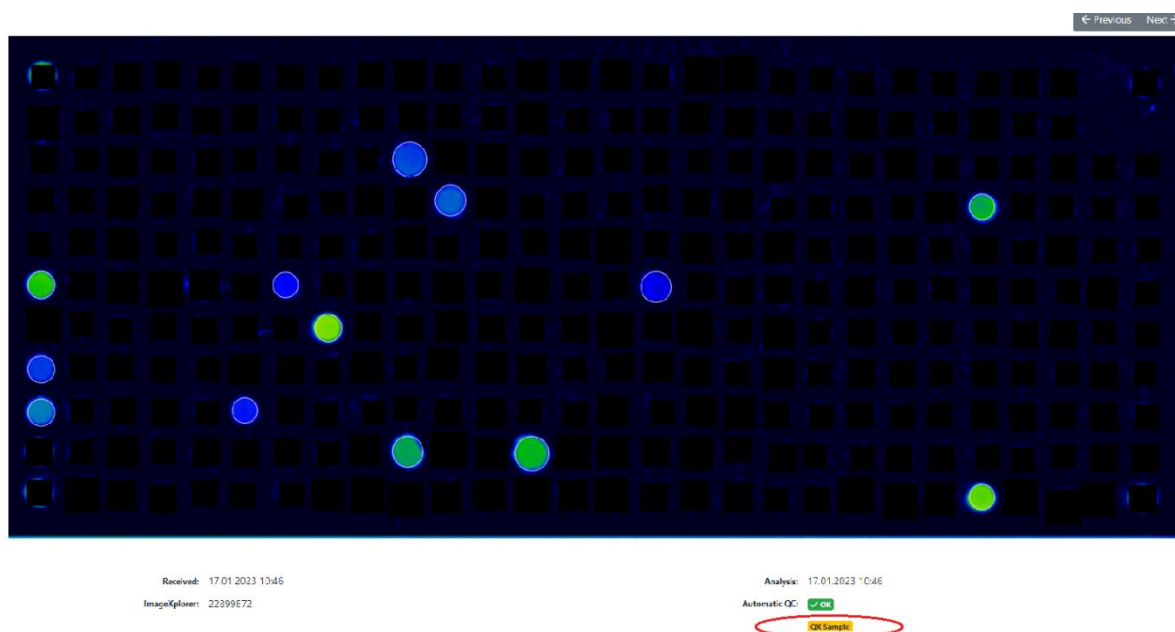
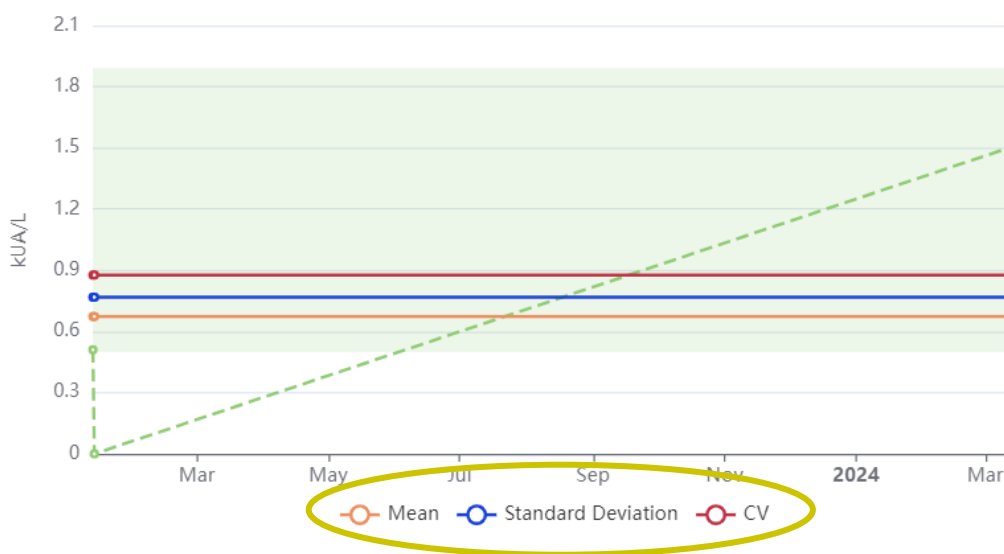
New Measurements

Search by barcode, patient, sample, tag, analysis date, a

Selected measurements ▾

- Download CSV format
- Download Excel format
- Approve
- Update interpretation
- Re-Process
- » Move
- Delete**

Tags	Product type
	FOX
	FOX
	ALEX2
	FOX
	ALEX2
	ALEX1

Delete measurements

You are about to delete the following measurements. Please note that this action is irreversible and cannot be undone.

Barcode	Analysis Date
02AAB160	16.12.2019 13:41

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Plant Food Food	M Jug r 3	0.40
Plant Food Food	M Jug r 1	0.38

Plant Food Food	M Jug r 3	0.40
Plant Food Food	M Jug r 1	0.38

Delete request scheduled now. Cancel delete request

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Sign in. Can be

Background jobs

View details about running background jobs and manually schedule runs.

[View dashboard](#)

Measurement delete / unapproval log & requests

Show logging and pending requests for measurement delete and unapproval.

[Show log & requests](#)

Measurement unapproval / delete requests

Requested UTC	Requesting User (Email)	QR-Code	Request Type	Action
11.10.2023 13:05:28	Test Test (Test@test.com)	02AYQ145	Unapprove	Accept Reject
11.10.2023 12:59:27	Test Test (Test@test.com)	02AYQ14C	Delete	Accept Reject

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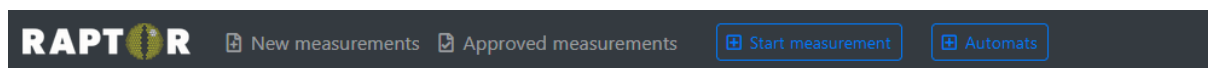
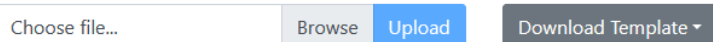
Add new user on Result Sharing Portal

General configuration

First name:

E-Mail address:

Username:

It is a good idea to use the users e-mail-address as username as well.

Status:

- Locked (user cannot log in)
- Enabled (user can log in)
- Locked (user cannot log in)

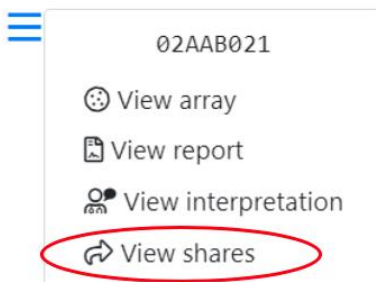
Last name:

Password:

Doctor Info Text:

This information will be printed on a patients report.

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≡ Share

You are about to share a measurement with a user of the User Portal. Please note that the doctor will have access to the results of the measurement. Once you share a measurement you'll be provided with a link to access the shared information via the encrypted User F

Barcode: 02AAB021

Share with: X Dr. (testuser2@macroarraydx.com)

If you can't find a specific user of the User Portal, go to the [tenant administration](#) to configure the list of available users of the User Portal.

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
The screenshot shows the RAPTOR interface with a dark header. Below the header, there are navigation tabs: 'New measurements', 'Approved measurements', 'Start measurement with ImageExplorer', and 'Automats'. The user is logged in as 'MADx Testtenant'. The main content area is titled 'Share' and contains a 'Share this measurement' button. Below this is a table with columns 'User', 'Link', and 'Shared'. A single row is visible with the user 'RP Bak', a long URL, and '7 months ago'. A red 'Revoke' button is located at the end of this row, with a yellow arrow pointing to it.

User	Link	Shared
RP Bak	https://resultshare.raptor-server.com/s/07PXEpC7ZNYthM4PdRfmct4NR5pUtWq5FJcqOY7Ry6oEx08USj	7 months ago

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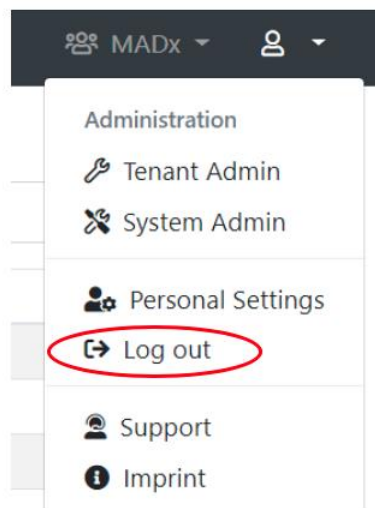Result Sharing Portal English (en) ▾

Welcome to the Result Sharing Portal

The Raptor-Server Result Sharing Portal allows users to view shared measurements. You will need a username and password to log in to view all measurements that have been shared with you. Also, a sharing link will allow to view specific shared measurements.

Please enter your username and your password:

Did you forget your password? No problem, you can [request a new one](#).

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Result Sharing Portal
English (en) ▼
Change password
Log out

Measurements

These are the measurements that are available for you. You may view the details of each measurement, or download the report PDF file.

Measurement	Product	Patient	Sample	Date
01AAK239	ALEX1		11852162	7/12/2019
01AAK281	ALEX1		04824821	12/18/2019
01AAK891	ALEX1		66560260	1/28/2020
02AAA015	ALEX2			2/21/2020

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Measurement 02AYQ143

Product: ALEX2
Patient:
Date of Birth:
Sample:
Info:
Doctor Info:
Sampling Date:
Assay Date: 10/12/2023
Processed Date: 10/12/2023
Approval Date: 10/12/2023

Report Language:

English (United Kingdom) ▾

You can request a new RAVEN interpretation that will be included in the report.

[Request new RAVEN interpretation](#)

You can download a PDF report that includes all information and results of this measurement.

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© Copyright by MacroArray Diagnostics