Q-Line **Grid**ION

Quick start guide

Q-Line

Pre-install

This Quick start guide sets up your Q-Line GridION™ for use.


See the following:

Q-Line GridION user manual
community.nanoporetech.com/to/gridion-q

Safety and regulatory information
community.nanoporetech.com/to/safety-q

For troubleshooting, see the user manual.

Q-Line GridION

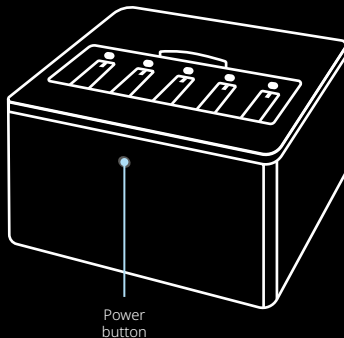**What's in
the box:**

Q-Line
GridION

White
Configuration
Test Cells (x5)

Power
cables*


DisplayPort to
HDMI adapter

* The Q-Line GridION ships with 5 x power cables (1 US, 1 UK, 1 EU, 1 CN, 1 AUS) for international use.

1

Set up your device

1. **Unpack** your Q-Line GridION device.
2. **Place the device** on a well-supported, strong, clean bench. Allow 30 cm clearance on all sides and top. **Do not cover ventilation grilles.** See user manual for detailed installation advice.
3. Attach the **cables** and **peripherals**.
4. Press the **power button** (shown on page 3).

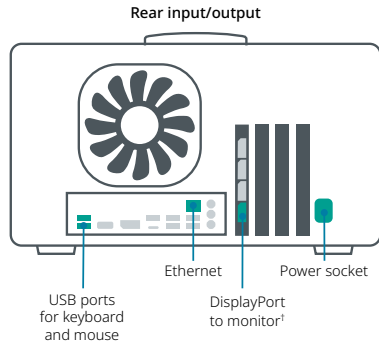¹If using HDMI only monitor, use the included DisplayPort to HDMI adapter.

Only the ports and connections shaded in green should be used.

2


Log in

1. To decrypt the **Master Key**, enter the default passphrase: **nanopore**
2. **Log in** to your Q-Line GridION
Password: **grid**
3. Update your Q-Line GridION **password**
Set a new password using the criteria below:
 - Minimum 8 characters
 - Must contain a capital letter and a lower case letter
 - Must contain a number or special character
 - Cannot be a palindrome (e.g. "radar")
 - Must differ from a previous password by more than 2 characters
4. Open the **Sequencing Software**
Click the wheel icon in the operating system favourites bar to load the Sequencing Software.

3

Perform a Position check

1. Slide the Q-Line GridION lid back, insert the CTCs under each clip in positions X1-X5 (as shown) and slide the lid shut.
2. Click **System** and then **Device information** in the Sequencing Software.
3. Click **Run Position check** and then **Start check** in the Sequencing Software.
4. Wait for the Sequencing Software to display  and '**Check complete. No issues detected**' to confirm the Position check has passed on your five flow cell positions*.
5. Remove your CTCs once your position check is complete.

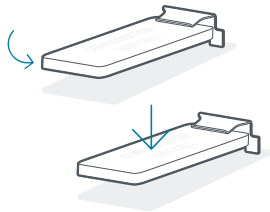*If any flow cell positions fail the Position check, ensure your CTCs are fully inserted into the flow cell positions, then repeat the Position check. Use the details in **Additional information** in this guide to contact Support if your Position check fails for any flow cell position **a second time**.

4

Power off the device

1. Click the menu under the power icon in the top right corner of the Ubuntu desktop for the drop-down menu.
2. Click **Power off** to power off your Q-Line GridION.

Note: Power off the Q-Line GridION at least once a month to ensure performance is not affected.

5

Access Q-Line documentation

Your Q-Line product information, resources, and documentation may be accessed at: nanoporetech.com/documentation.

Legacy content requires a user log-in

Additional information

View Store

Oxford Nanopore Q-Line products can be found here: store.nanoporetech.com

Note: Legacy content requires a user log-in.

Software updates

Oxford Nanopore will inform you of Sequencing Software updates.

DO NOT update your Q-Line GridION with MinKNOW from the Nanopore Community.

Documentation

Documentation may be accessed at nanoporetech.com/documentation

Note: Legacy content requires a user log-in.

Support

Technical Support will contact you shortly to ensure the device is set up correctly.

Please contact support@nanoporetech.com if you have any issues during setup.

Visit nanoporetech.com/support for information, live chat, and phone assistance.

Warranty

Extend the license and warranty for your device here:

nanoporetech.com/products/services/oxford-nanopore-support-plans

For flow cell warranty, visit: community.nanoporetech.com/to/warranty

Recycle used flow cells

Oxford Nanopore is committed to environmental sustainability. You can help by sending your flow cells for recycling. Find out how: community.nanoporetech.com/support/returns

Technical specification

Model number	Q-GRD-MK1	
Supply voltage (V)	100-240 AC $\pm$ 10% (50-60 Hz)	
Maximum rated current (A)	6.5	
Maximum rated power (W)	650	
Size (H x W x D) (mm)	220 x 365 x 370	
Weight (kg)	14.4	
Installation ports	1 x Ethernet port (1 Gbps) 1 x HDMI/DisplayPort to monitor 1 x USB-A for keyboard	1 x USB-A for mouse 1 x Power cable
Software installed	Ubuntu, GridION OS, Sequencing Software	
Compute specification	7 TB SSD storage, 64 GB RAM, minimum 8 core Intel CPU, 1 x NVIDIA GPU	
Environmental conditions	Functional range of electronics is within environmental temperatures of +5°C to +40°C. Users should allow 30 cm clearance to the rear and sides of the device. Designed to sequence in environmental temperatures of +18°C to +25°C. Intended for indoor use. Can be used up to altitudes of 2,000 m. Use within 30%–75% relative non-condensing humidity limits. The device has a Pollution Degree 2. This is an Overvoltage 2 device. WARNING: Rear of instrument heats up during operation. WARNING: Use the mains power cable supplied, or a suitable, approved, and rated power cable. Ensure the device is connected to a protective earth mains supply.	

Oxford Nanopore Technologies

phone +44 (0)845 034 7900

email support@nanoporetech.com

 [oxford-nanopore-technologies](https://www.linkedin.com/company/oxford-nanopore-technologies)

 [@nanopore](https://twitter.com/nanopore)

 [@nanoporetech.com](https://www.facebook.com/nanoporetech.com)


Gosling Building,
Edmund Halley Road,
Oxford Science Park,
Oxfordshire,
OX4 4DQ, UK

www.nanoporetech.com

Oxford Nanopore Technologies, the Wheel icon, GridION and MinKNOW, are registered trademarks or the subject of trademark applications of Oxford Nanopore Technologies plc in various countries. Information contained herein may be protected by copyright, patents or patents pending of Oxford Nanopore Technologies plc. All other brands and names contained are the property of their respective owners. © 2026 Oxford Nanopore Technologies plc. All rights reserved. Oxford Nanopore Technologies products are RUO. Products labelled/branded as Oxford Nanopore Diagnostics may be RUO or may be regulated as in-vitro diagnostic devices in some jurisdictions, please check individual product labelling.

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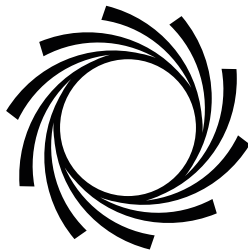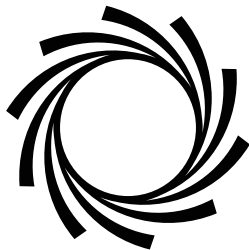
community.nanoporetech.com/to/gridion-q

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community.nanoporetech.com/to/safety-q

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